

Description	Deep-Well Microplates			
Biotix Part #	DP-1200-9CV	DP-1200-9CVS	DP-2200-9CV	DP-2200-9CVS
Product Description	96-square well	96-square well	96-square well	96-square well
Maximum Volume	1.2 ml	1.2 ml	2.2 ml	2.2 ml
Certified Pre-Sterile	No	Yes	No	Yes
Composition	Medical-grade virgin polypropylene			
Features	Plate dimensions conform to ANSI/SBS 1-2004 standards for automated workstations and high throughput processing			
Packaging	5 plates/pack, 10 packs/case			
Physical Data	DNase, RNase and Endotoxin (pyrogen) free			





1.2 ml



2.2 ml

## Product Dimensions

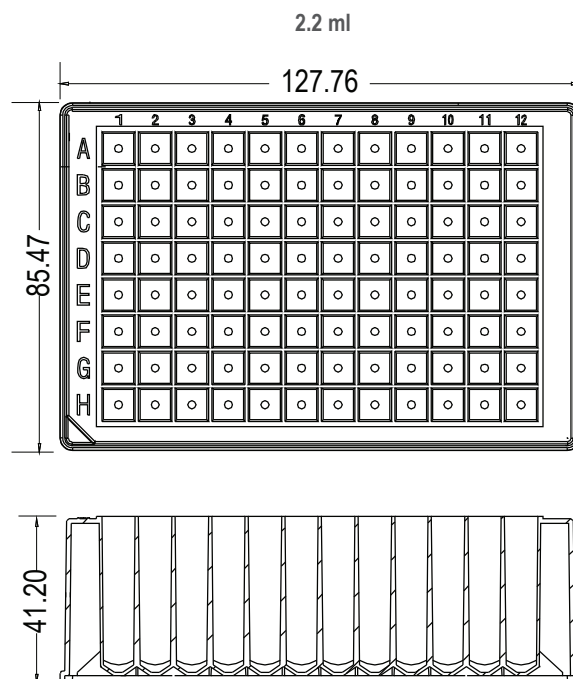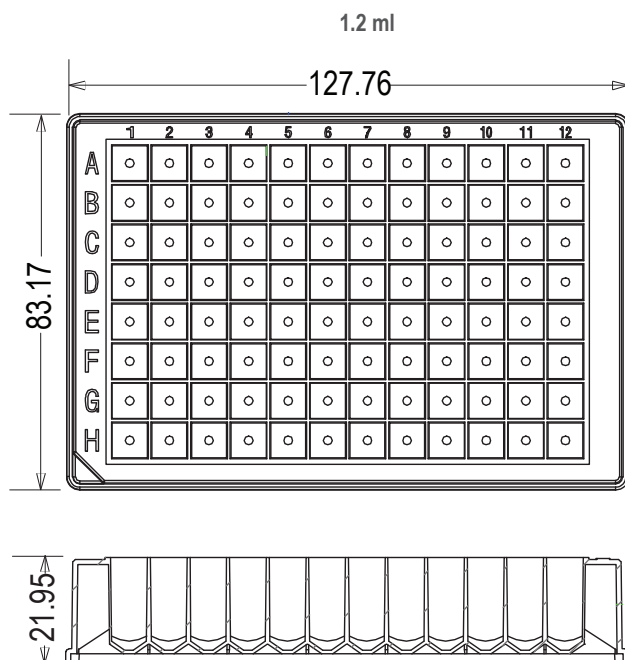

## Quality Testing

<b>RNase/DNase</b>	Products are washed in distilled water and concentrated via centrifugation. Samples are added to previously established nucleic acid standards, incubated for one hour at 37°C, and tested on a 2% gel using electrophoresis. Products must show no degradation of standards to pass. Test sensitivity is 10 <sup>-7</sup> Kunitz units/μl.
<b>Nucleic Acid</b>	Products are washed in distilled water and concentrated via centrifugation. Then, samples are added to protocol specified PCR reactions and thermal cycled for 50 cycles. A 2% agarose gel electrophoresis is used to examine experimental samples, positive controls, and negative controls. To pass, product samples must show no DNA amplification. Test sensitivity is 10 ng.
<b>Endotoxin/Pyrogen</b>	Products are tested for endotoxins by using the Limulus Amebocyte Lysate (LAL) gel assay according to FDA guidelines. Test sensitivity is 0.06 EU/ml.
<b>Trace Metal</b>	Products are washed in distilled water. The sample is then tested using reflectometry using a single strip test for the following metals: Ca, Cu, Fe, K, Mg, Mn and Ni. Standard solutions are used as positive controls. A reader is used to detect metals to a sensitivity of 500 mg/L.
<b>PCR Inhibitor</b>	Products are tested via PCR amplification and gel electrophoresis analysis. Samples must show normal amplification to be considered free of PCR inhibitors.
<b>Sterilization</b>	Products are sterilized using electron beam irradiation.

©2013, Biotix, Inc. All rights reserved

Biotix, Fluid Innovation, Delta Filter, X-Resin, Blade, FlexFit, Statix, Leave Nothing Behind, BioReady, Tip Eject, and XTreme Quality are registered trademarks or trademarks of Biotix, Inc. All other brands and names contained herein are the property of their respective owners.