


xTIP4 Tip Eject™ Reload

Experience convenience and quality
while reducing your lab's waste by 30%

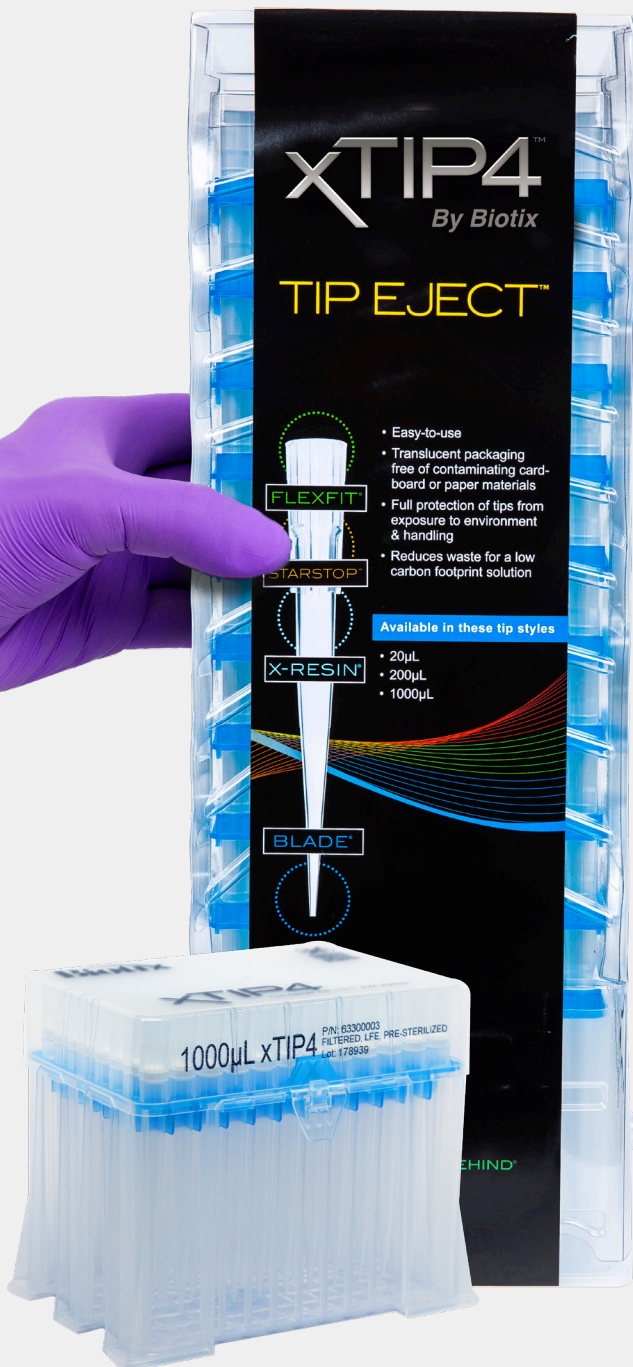


This self-aligning reload solution allows you to easily reload your racks without coming into contact with the tips. The rigid plastic covering also shields unused tips from contamination for later use.

- Available only for LTS and xPIPETTE compatible tips.
- Vertically stacked snapcards take up less space on your lab bench.
- Reduces plastic waste by 30% when compared to racked tips.

Pipette tip racks are fully autoclaveable at 121° C after transfer from reload.



Learn more about our sustainability initiatives at BIOTIX.com

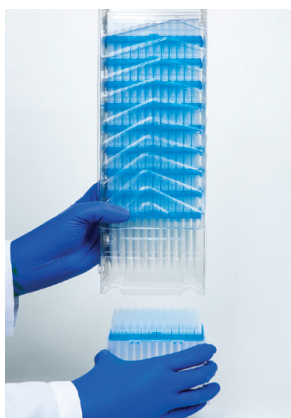

Step 1: Remove the base

Remove the lower protective rack base.



Step 2: Align empty rack

Align and place Tip Eject over your empty Biotix rack.



Step 3: Push down

Push down until a "click" is heard, indicating tips are locked into rack.



Step 4: Replace stack over empty rack base

Repeat to refill empty racks as needed

Catalog No.	Description	Packaging
-------------	-------------	-----------

xTIP4 Tip Eject Reload, Compatible with xPIPETTE and LTS pipettes

63300015	0.2-20 μ L Sterilized	10 refills/pack
63300016	20-200 μ L Sterilized, Low Retention	
63300018	100-1000 μ L Sterilized, Low Retention	8 refills/pack
63300019	100-1200 μ L Sterilized, Low Retention	

xTIP4 Empty Racks

63300151	Empty Racks for Tip Eject Reloads, 20-300 μ L tips	10 racks/pack
63300152	Empty Racks for Tip Eject Reloads, 1000-1200 μ L tips	

