

# Biotix Tip Eject Reload

Compatible with Biotix xTIP® and Rainin LTS™ tips

Experience convenience and quality while reducing your lab's waste by 30%




This self-aligning reload solution allows you to easily reload your racks without coming into contact with the tips. The rigid plastic covering also shields unused tips from contamination for later use.

- Available only for LTS and xPIPETTE compatible tips.
- Vertically stacked snapcards take up less space on your lab bench.
- Reduces plastic waste by 30% when compared to racked tips.

Pipette tip racks are fully autoclaveable at 121° C after transfer from reload.



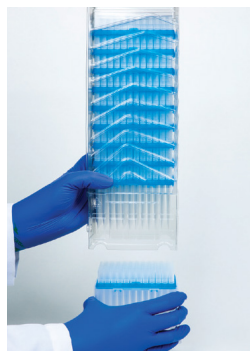

**Step 1: Remove the base**

Remove the lower protective rack base.



**Step 2: Align empty rack**

Align and place Tip Eject over your empty Biotix rack.



**Step 3: Push down**

Push down until a “click” is heard, indicating tips are locked into rack.



**Step 4: Replace stack over empty rack base**

Repeat to refill empty racks as needed

Volume $\mu$ L	Pre-Sterilized	Tip Eject Reload	Tip Eject Packaging
10	Yes		
	No		
10 XL	Yes		
	No		
20	Yes	76266-128	10 refills/pack
	No		
200	Yes		
	No		
250	Yes	76266-130	10 refills/pack
	No		
300	Yes		
	No		
1000	Yes	76266-134	8 refills/pack
	No		
1200	Yes	76266-136	8 refills/pack
	No		