

Experience convenience and quality while reducing your lab's waste by 30%



This self-aligning reload solution allows you to easily reload your racks without coming into contact with the tips. The rigid 100% recycled plastic covering shields unused tips from contamination for later use.

- Available for LTS and xPIPETTE compatible tips.
- 100% recycled packaging material
- Easy-to-use fast reload system
- Reduces plastic waste by 30% when compared to racked tips.
- Vertically stacked snapcards take up less space on your lab bench.

Pipette tip racks are fully autoclaveable at 121° C after transfer from reload.



Learn more about our sustainability initiatives at [BIOTIX.com](https://www.biotix.com)



**Step 1: Remove the base**

Remove the lower protective rack base.



**Step 2: Align empty rack**

Align and place Tip Eject over your empty Biotix rack.



**Step 3: Push down**

Push down until a “click” is heard, indicating tips are locked into rack.



**Step 4: Replace stack over empty rack base**

Repeat to refill empty racks as needed

Volume $\mu\text{L}$	Sterilized	xTIP Tip Eject Reload	xTIP Tip Eject Packaging
20	Yes	63300015	10 refills/pack
250	Yes	63300016	10 refills/pack
300	Yes	63305951	10 refills/pack
1000	Yes	63300018	8 refills/pack
1200	Yes	63300019	8 refills/pack

Volume $\mu\text{L}$	Empty Rack	Packaging
10-300	63300151	10 Rack/Pack
1000-1200	63300152	10 Rack/Pack